



## *General Protocol & Storage*

### Product Description

A set of Stellaris RNA FISH Probes is comprised of up to 48 singly labeled oligonucleotides designed to selectively bind to targeted transcripts. Stellaris RNA FISH Probes bound to target RNA produce fluorescent signals that permit detection of single RNA molecules as diffraction-limited spots by conventional fluorescence microscopy.

### Storage Guidelines

#### *Stellaris RNA FISH Probes*

Stellaris RNA FISH Probes are shipped dry and can be stored at +2 to +8 °C in this state. Dissolved probe mix should be subjected to a minimum number of freeze-thaw cycles. For daily and short-term use of dissolved probe mix, storage at +2 to +8 °C in the dark for up to a month is recommended. For storage lasting longer than a month, we recommend aliquoting and freezing probes in the dark at -15 to -30 °C.

#### *Stellaris RNA FISH Hybridization Buffer*

Stellaris RNA FISH Hybridization Buffer should be stored at +2 to +8 °C for short-term and long-term use.

#### *Stellaris RNA FISH Wash Buffer A and Wash Buffer B*

Stellaris RNA FISH Wash Buffers A and B should be stored at room temperature for short-term and long-term use.

### Reagents and Equipment

#### *Reagents and Consumables:*

- a) TE buffer (10 mM Tris-HCl, 1 mM EDTA, pH 8.0)
- b) 20% Electron Microscopy Grade Formaldehyde
- c) 10X Phosphate Buffered Saline (PBS), RNase-free
- d) Nuclease-free water
- e) Ethanol for molecular biology
- f) Triethanolamine
- g) Acetic Anhydride
- h) Chloroform
- i) Deionized Formamide
- j) Stellaris RNA FISH Hybridization Buffer (LGC Biosearch Technologies Cat# SMF-HB1-10)
- k) Stellaris RNA FISH Wash Buffer A (LGC Biosearch Technologies Cat# SMF-WA1-60)
- l) Stellaris RNA FISH Wash Buffer B (LGC Biosearch Technologies Cat# SMF-WB1-20)
- m) 4',6-diamidino-2-phenylindole (DAPI)
- n) Prolong<sup>®</sup> Gold Antifade Mountant (ThermoFisher Scientific Cat #P36930)
- o) 24 x 60 mm rectangular coverglass
- p) RNase free consumables such as pipette tips
- q) Humidified chamber (or equivalent): 150 mm tissue culture plate; a single water-saturated paper towel placed alongside the inner chamber edge
- r) 37 °C laboratory oven

### Microscope:

- Wide-field fluorescence microscope (e.g., Nikon Eclipse Ti or equivalent). We provide limited support for confocal applications.
- A high numerical aperture (>1.3) and 60-100x oil-immersion objective.
- Strong light source, such as a mercury or metal-halide lamp (newer LED-based light sources may also be sufficient).
- Filter sets appropriate for the fluorophores.
- Standard cooled CCD camera, ideally optimized for low-light level imaging rather than speed (13 µm pixel size or less is ideal).

### Preparation of Reagents

**NOTE:** When performing Stellaris RNA FISH, it is imperative to limit RNA degradation. Please ensure that all consumables and reagents are RNase-free. Recipes below are for set volumes. Please adjust accordingly.

#### Reconstituting the dried probe stock:

**ShipReady Probe Set (1 nmol):** A ShipReady probe set can provide up to 80 hybridizations. Re-dissolve the dried oligonucleotide probe blend in 80 µL of TE buffer (10 mM Tris-HCl, 1 mM EDTA, pH 8.0) to create a probe stock of 12.5 µM. *Mix well by pipetting up and down*, and then vortex and centrifuge briefly.

**DesignReady or Custom Probe Set (5 nmol):** A DesignReady or custom probe set can provide up to 400 hybridizations. Re-dissolve the dried oligonucleotide probe blend in 400 µL of TE buffer (10 mM Tris-HCl, 1 mM EDTA, pH 8.0) to create a probe stock of 12.5 µM. *Mix well by pipetting up and down*, and then vortex and centrifuge briefly.

#### Fixation Buffer:

*Final composition is 3.7% (vol./vol.) formaldehyde in 1X PBS*

**For a final volume of 50 mL, mix:**

- 1 mL 20% Formaldehyde solution
- 5 mL 10X Phosphate Buffered Saline (PBS), RNase-free
- 35 mL Nuclease-free water

#### TEA Buffer (10X):

**NOTE:** *This buffer is light sensitive. It is suggested to wrap reagent bottle in foil.*

**For a final volume of 100 mL, mix:**

- 13.3 mL Triethanolamine
- 60 mL Nuclease-free water H<sub>2</sub>O
- Then bring to pH 8.0 using HCL (~5-10 mL of HCL)
- Once at pH 8.0, add water to 100 mL final volume

#### Hybridization Buffer:

*Final composition is 10% (vol./vol.) formamide in Hybridization Buffer*

**Hybridization Buffer should be mixed fresh for each experiment:**

Due to viscosity of the solution, we recommend accounting for a 10% final volume excess in order to have enough Hybridization Buffer for all of your samples.

**For a final volume of 1 mL, mix:**

- 900 µL Stellaris RNA FISH Hybridization Buffer (LGC Biosearch Technologies Cat# SMF-HB1-10)
- 100 µL Deionized Formamide

**NOTE:** Do not freeze Hybridization Buffer.

**WARNING!** Formamide is a teratogen that is easily absorbed through the skin and should be used in a chemical fume hood. Please consult the appropriate SDS (Safety Data Sheet) prior to use.

**WARNING!** Be sure to let the formamide warm to room temperature before opening the bottle.

**WARNING!** Formaldehyde is a known human carcinogen and should be use in a chemical fume hood. Please consult the appropriate SDS (Safety Data Sheet) prior to use.

### **Wash Buffer A (50 mL):**

Final composition is 10% (vol./vol.) formamide in 1X Wash Buffer A

#### **Mix and dilute Wash Buffer A fresh for each experiment:**

##### **For a final volume of 50 mL, mix:**

- 10 mL Stellaris RNA FISH 5X Wash Buffer A (LGC Bioscience Technologies Cat# SMF-WA1-60)
- Add 5 mL Deionized Formamide
- Add 35 mL Nuclease-free water
- Mix well by vortexing gently.

### **Wash Buffer B:**

#### **Add Nuclease-free water to Wash Buffer B bottle upon first use.**

Add 88 mL of Nuclease-free water to bottle (LGC Bioscience Technologies Cat# SMF-WB1-20) before use. Mix thoroughly.

### **Nuclear Stain for use after hybridization:**

4',6-diamidino-2-phenylindole (DAPI) prepared in Wash Buffer A (see above) at 5 ng/mL. This solution is to be used in Step J below.

### **Mounting media:**

Prolong Gold Antifade Mountant (ThermoFisher Scientific Cat #P36930).

**NOTE:** Samples mounted with Prolong Gold should be allowed to cure overnight at room temperature and then imaged the following day.

## ***Protocol for Fresh Frozen Mouse Brain Tissue***

### **Day 1- Slide Prep and Hybridization**

1. Thaw the slide-mounted tissue section to room temperature.
2. Immerse the slide in cold 4% E.M. grade paraformaldehyde in 1X PBS for 15 minutes.
3. Wash twice with 1X PBS for 5 minutes.
4. Dip the slide in nuclease-free water.
5. Dip the slide in 1X TEA buffer.
6. Immerse the slide in 1X TEA + Acetic Anhydride for 10 minutes (Stirring!) \*\*\*

**Note:** Acetic Anhydride must be added fresh each time. Add Acetic Anhydride to fresh 1X TEA buffer after the second 1X PBS wash in step 3, about 1 minute before dipping slides in DEPC water. For 50 mL of 1X TEA, add 63  $\mu$ L of Acetic Anhydride.

7. Immerse the slide in 2X SSC for 3 minutes.
8. Immerse the slide in 70% ethanol for 3 minutes.
9. Immerse the slide in 95% ethanol for 3 minutes.
10. Immerse the slide in 100% ethanol for 3 minutes.
11. Immerse the slide in Chloroform for 5 minutes.
12. Immerse the slide in 100% ethanol for 3 minutes.
13. Immerse the slide in 95% ethanol for 3 minutes.
14. Let air dry for 90+ minutes (but no longer than 4 hours).

### **Hybridization in frozen tissue sections**

If frozen before using, warm the reconstituted probe solution to room temperature. Mix well by vortexing, then centrifuge briefly.

To prepare the hybridization solution, add 4.0  $\mu$ L of probe stock solution to 200  $\mu$ L of hybridization buffer, and then vortex and centrifuge. This creates a working probe solution of 250 nM. This solution will be used on step 16.

15. Assemble a humidified chamber: 150 mm tissue culture plate; a single water-saturated paper towel placed alongside the inner chamber edge. This chamber will help prevent evaporation of the probe solution from the tissue section.
16. After slide has dried for 90+ minutes, dispense 200  $\mu$ L of hybridization buffer containing probe onto the tissue sections of the slide.

17. Carefully place a clean 24 x 60 mm rectangular coverglass over the hybridization solution to completely cover the tissue sections and allow for even distribution of the hybridization solution. Place the slide in the humidified chamber, cover with the tissue culture lid, and seal chamber with parafilm.
18. Incubate in the dark at 37 °C for at least 4 hours (incubation can be continued up to 16 hours).

## Day 2- In situ washes

19. Immerse the slide in wash buffer A, and allow the submerged coverglass to slide off the tissue section. Gentle agitation may be required to remove the coverglass.
20. Incubate in the dark at 37 °C for 30 minutes.
21. Decant wash buffer A, and then add DAPI nuclear stain (wash buffer consisting of 5 ng/mL DAPI) to counterstain the nuclei.
22. Incubate in the dark at 37 °C for 30 minutes.
23. Decant DAPI staining buffer, and then immerse slide in Wash Buffer B for 3 minutes.
24. Immerse slide in 50% ethanol for 3 minutes.
25. Immerse slide in 85% ethanol for 3 minutes.
26. Immerse slide in 100% ethanol for 3 minutes.
27. Let air dry for 5-10 minutes.
28. Add a drop or two (approximately 50-100 µL) of Prolong Gold Antifade Mountant onto the tissue sections. Cover with a clean 24 x 60 mm coverglass, allowing the antifade to spread evenly across the tissue sections.
29. Allow Prolong Gold to cure overnight, in the dark, at room temperature.
30. Seal the coverglass perimeter with clear nail polish, and allow to dry in the dark.
31. Proceed to imaging.

## References

1. Raj, A., van den Bogaard, P., Rifkin S.A., van Oudenaarden, A., and Tyagi, S. Imaging individual mRNA molecules using multiple singly labeled probes. *Nat. Methods*. 2008; 5: 877-9
2. Femino, A.M., Fay, F.S., Fogarty, K., and Singer, R.H. Visualization of single RNA transcripts in situ. *Science* 1998; 280: 585-90
3. Raj A, Tyagi S. Detection of individual endogenous RNA transcripts in situ using multiple singly labeled probes. *Methods in Enzymology* 2010; 472: 365-86.

## Guidelines for Citing the Use of Stellaris RNA FISH Probes and Methods in Scientific Publications

Please acknowledge the use of Stellaris RNA FISH Probes and/or protocols in the experimental **Materials and Methods** or **Methods** section of your manuscript. Refer to the following examples as guidelines for proper citation of the Stellaris RNA FISH probe sets and/or protocols:

**Citing ShipReady and DesignReady Probe Sets:** "Stellaris® FISH Probes recognizing <catalogued gene set name> and labeled with <your dye of choice> (Catalog #, Biosearch Technologies, Inc., Petaluma, CA) were hybridized to <samples>, following the manufacturer's instructions available online at [www.biosearchtech.com/stellarisprotocols](http://www.biosearchtech.com/stellarisprotocols). Briefly, <describe any deviations from the published protocol or a short summary of what was actually performed>."

**Citing Custom Probe Sets Designed with the Stellaris® FISH Probe Designer:** "Custom Stellaris® FISH Probes were designed against <your RNA of interest (include NM# and nucleotides covered if relevant)> by utilizing the Stellaris® FISH Probe Designer (Biosearch Technologies, Inc., Petaluma, CA) available online at [www.biosearchtech.com/stellarisdesigner](http://www.biosearchtech.com/stellarisdesigner). The <samples> were hybridized with the <your RNA of interest> Stellaris FISH Probe set labeled with <your dye of choice> (Biosearch Technologies, Inc.), following the manufacturer's instructions available online at [www.biosearchtech.com/stellarisprotocols](http://www.biosearchtech.com/stellarisprotocols). Briefly, <describe any deviations from the published protocol or a short summary of what was actually performed>."

**Citing Custom Probe Sets Utilizing Previously Published Sequences:** "Custom Stellaris® FISH Probes recognizing <your RNA of interest (include NM# and nucleotides covered if relevant)> and labeled with <your dye of choice>, were purchased from Biosearch Technologies, Inc. (Petaluma, CA). Probe set sequences utilized in the experiments have been previously described <cite published manuscript>. The <samples> were hybridized with the <your RNA of interest> Stellaris® FISH Probe set, following the manufacturer's instructions available online at [www.biosearchtech.com/stellarisprotocols](http://www.biosearchtech.com/stellarisprotocols). Briefly, <describe any deviations from the published protocol or a short summary of what was actually performed>."

**Citing 3' Amine Oligos in Plates Used for Stellaris RNA FISH Designed with the Stellaris® FISH Probe Designer:** "Custom 3' amine oligos in plates were designed against <your RNA of interest (include NM# and nucleotides covered if relevant)> by utilizing the Stellaris® FISH Probe Designer (Biosearch Technologies, Inc., Petaluma, CA) available online at [www.biosearchtech.com/stellarisdesigner](http://www.biosearchtech.com/stellarisdesigner). Probes were labeled with <your dye of choice> using <insert your labeling protocol or citation of previously published labeling protocol>. The <samples> were hybridized with the <your RNA of interest> oligonucleotides (Biosearch Technologies, Inc.), following the manufacturer's instructions available online at [www.biosearchtech.com/stellarisprotocols](http://www.biosearchtech.com/stellarisprotocols). Briefly, <describe any deviations from the published protocol or a short summary of what was actually performed>."

**Citing 3' Amine Oligos in Plates Used for Stellaris RNA FISH Using Previously Published Sequences:** "Custom 3' amine oligos in plates recognizing <your RNA of interest (include NM# and nucleotides covered if relevant)> were purchased from Biosearch Technologies, Inc. (Petaluma, CA). Probe set sequences utilized in the experiments have been previously described <cite published manuscript>. Probes were labeled with <your dye of choice> using <insert your labeling protocol or citation of previously published labeling protocol>. The <samples> were hybridized with the <your RNA of interest> oligonucleotides (Biosearch Technologies, Inc.), following the manufacturer's instructions available online at [www.biosearchtech.com/stellarisprotocols](http://www.biosearchtech.com/stellarisprotocols). Briefly, <describe any deviations from the published protocol or a short summary of what was actually performed>."

## Technical Support

If you require additional information or technical assistance please feel free to e-mail our Technical Support Group at: [techsupport@biosearchtech.com](mailto:techsupport@biosearchtech.com). Our knowledgeable staff is also available for telephone consultation from 8:00 AM to 5:00 PM, Monday through Friday, Pacific Time. Please contact us at:

1.800.GENOME.1 (436.6631) US & Canada only

+1.415.883.8400 telephone

+1.415.883.8488 fax

2199 S. McDowell Blvd, Petaluma, CA 94954-6904 USA | [info@biosearchtech.com](mailto:info@biosearchtech.com)

©2016 Biosearch Technologies, Inc. All rights reserved.

This product is sold under license from Rutgers, The State University of New Jersey, and may be used under its patent rights for Research Use Only. Information on licensing programs for the commercial and/or diagnostic use of these products is available by contacting [licensing@biosearchtech.com](mailto:licensing@biosearchtech.com).

Biosearch, the Biosearch Technologies logo, and Stellaris are registered trademark of Biosearch Technologies, Inc., Products and technologies appearing in this brochure may have trademark or patent restrictions associated with them. Please see [www.biosearchtech.com/legal](http://www.biosearchtech.com/legal) for full legal disclosure.

UI-207727 Rev. 1.0  
Eff. Date: 03 Feb 2016