## BIOSEARCH TECHNOLOGIES Advancing Nucleic Acid Technology<sup>SM</sup>

## (FISH) ASSAY

- and in tissue.



with clear nail polish.



# Improving Dye Brightness and Photostability in Stellaris<sup>®</sup> RNA FISH Assays

Mary Katherine Johansson, Sally R. Coassin, Arturo V. Orjalo, Matt H. Lyttle and Ron M. Cook 81 Digital Drive, Novato, CA 94949-5728 USA

Next, a GAPDH Stellaris probe set was prepared with the nitrobenzyl-Quasar 670 adduct. Unfortunately, compared to a standard GAPDH Quasar 670 probe set, consistent spots of mRNA were not seen throughout the slide. This suggests that the TSQ moiety is too hydrophobic and interfered with target hybridization.

### CONCLUSIONS

- Both diffusional antifades and covalent TSQ can improve dye brightness and photostability.
- The covalent TSQ moieties worked well in test poly-T sequences but appeared to interfere with hybridization in a Stellaris assay with a mixed base probe set.
- Efforts are ongoing to find an antifade as effective and reproducible as Vectashield.

Glucose oxidase catalyzes:  $\beta$ -D-glucose + O<sub>2</sub>  $\rightarrow$  D-gluconolactone + H<sub>2</sub>O<sub>2</sub> \*\*\* GLOX is incompatible with pH dependent dyes such as fluorescein.

Adding GLOX to a solution with FAM makes the fluorescence disappear.

- Dye bleaching occurs within seconds with FAM and Quasar 670 dyes under standard wide field microscopy conditions using a metal halide lamp.
- Several potential triplet state quenchers and antioxidants were added to the final mounting buffer solution to test antifade efficacy.
- Dye fluorescence intensity over time (photostability) and initial fluorescence intensity (brightness) were calculated for both FAM and Quasar 670 dyes.
- The antifade efficacy depends on several parameters including concentration, pH, viscosity, and buffer composition.
- Vectashield<sup>®</sup> from Vector Laboratories is a very effective antifade for both FAM and Quasar 670 dyes.
- Some other antifade compositions were almost as effective as Vectashield, but not as reproducible.

- **2012**, *9*, 68-71.



#### REFERENCES

• M. Batish, A. Raj, S. Tyagi, Single molecule imaging of RNA in situ, Methods Mol. Biol., 2011, 714, 3-13. • R. B. Altman et al., Cyanine fluorophore derivatives with enhanced photostability, Nature Methods,

• M. Levitus, S. Ranjit, Cyanine dyes in biophysical research: the photophysics of polymethine fluorescent dyes in biomolecular enviornments, Quarterly Reviews of Biophysics, 2011, 44, 123-151.

itellaris®, CAL Fluor®, and Quasar® are registered trademarks of Biosearch Technologies. Inc

see www.biosearchtech.com/legal for full legal disclosu his product is sold under license from the University of Medicine and Dentistry of New Jersey and may be used under its patent rights for Research Use Only.

Scan this QR Code to learn more about Stellaris RNA FISH Probes.

