

# Product insert

## Charité 2019-nCoV BHQ ValuPanel Reagents

For Research Use Only. Not for use in diagnostic procedures.

### Products

Size	Cat no.	Product description	Delivered amount
Small	RdRP-SARSr-F-20	Charite RdRP_SARSr Forward Primer	20 nmol
	RdRP-SARSr-R-20	Charite RdRP_SARSr Reverse Primer	20 nmol
	RdRP-SARSr-BHQP1-5	Charite RdRP_SARSr Pan Sarbeco FAM-BHQ Probe	5 nmol
	RdRP-SARSr-BHQP2-5	Charite RdRP_SARSr 2019-nCoV FAM-BHQ Probe	5 nmol
	E-Sarbeco-F-20	Charite E_Sarbeco Forward Primer	20 nmol
	E-Sarbeco-R-20	Charite E_Sarbeco Reverse Primer	20 nmol
	E-Sarbeco-BHQP-5	Charite E_Sarbeco FAM-BHQ Probe	5 nmol
Medium	RdRP-SARSr-F-100	Charite RdRP_SARSr Forward Primer	100 nmol
	RdRP-SARSr-R-100	Charite RdRP_SARSr Reverse Primer	100 nmol
	RdRP-SARSr-BHQP1-25	Charite RdRP_SARSr Pan Sarbeco FAM-BHQ Probe	25 nmol
	RdRP-SARSr-BHQP2-25	Charite RdRP_SARSr 2019-nCoV FAM-BHQ Probe	25 nmol
	E-Sarbeco-F-100	Charite E_Sarbeco Forward Primer	100 nmol
	E-Sarbeco-R-100	Charite E_Sarbeco Reverse Primer	100 nmol
	E-Sarbeco-BHQP-25	Charite E_Sarbeco FAM-BHQ Probe	25 nmol
Large	RdRP-SARSr-F-1000	Charite RdRP_SARSr Forward Primer	1000 nmol
	RdRP-SARSr-R-1000	Charite RdRP_SARSr Reverse Primer	1000 nmol
	RdRP-SARSr-BHQP1-250	Charite RdRP_SARSr Pan Sarbeco FAM-BHQ Probe	250 nmol
	RdRP-SARSr-BHQP2-250	Charite RdRP_SARSr 2019-nCoV FAM-BHQ Probe	250 nmol
	E-Sarbeco-F-1000	Charite E_Sarbeco Forward Primer	1000 nmol
	E-Sarbeco-R-1000	Charite E_Sarbeco Reverse Primer	1000 nmol
	E-Sarbeco-BHQP-250	Charite E_Sarbeco FAM-BHQ Probe	250 nmol

The Charité 2019-nCoV BHQ ValuPanel Reagents include FAM-BHQ probes and primers delivered in individual tubes in dried format using sequences designated by the Charité protocol for 2019-nCoV. For Research Use Only. Not for use in diagnostic procedures.

We offer this set of ValuPanel Reagents in three different sizes, small, medium, and large, where customers can select which probes and primers are required for their needs. The number of reactions will depend on the protocol selected by the customer.

### Storage

Charité 2019-nCoV BHQ ValuPanel Reagents are shipped dry at ambient temperature. They may be stored at +2 to +8 °C in this state.

*Continue for more information»*



# Product insert

## Charité 2019-nCoV BHQ ValuPanel Reagents

For Research Use Only. Not for use in diagnostic procedures.

---

### Certificates of Analysis (CoA)

Please contact [techsupport@lgcgroup.com](mailto:techsupport@lgcgroup.com) to obtain a customer-specific CoA for each ordered probe and primer.

**Integrated tools.  
Accelerated science.**

   @LGCBiosearch | [biosearchtech.com](http://biosearchtech.com)

All trademarks and registered trademarks mentioned herein are the property of their respective owners. All other trademarks and registered trademarks are the property of LGC and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details. No part of this publication may be reproduced or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording or any retrieval system, without the written permission of the copyright holder. © LGC Limited, 2020. All rights reserved. GEN782/GL/0420

**BIOSEARCH™**  
**TECHNOLOGIES**  
GENOMIC ANALYSIS BY LGC